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Tips for Keeping your Laboratory Freezer Running Efficiently

Monitor Temperatures

Installing an alarmed thermometer monitoring system, like Heathrow Scientific's Freezer Thermometer, can alert you early to any fluctuations in temperatures that could indicate the need for cleaning or maintenance.

Organize to Ensure Proper Air Flow

Our Freezer Rack systems not only use all optimal storage space available; they also allow for air flow through the unit. With materials stacked and organized, it makes it easier to locate samples quickly, greatly reducing the amount of time the door remains open.

Remove Ice and Frost Routinely

An accumulation of ice and frost on racks and doors can interrupt air flow and overwork the freezer. Defrost on a routine basis to keep your unit in proper working order.

Clear Debris from Compressor Coils

Vacuum or use a brush to clean the compressor coils on the outside of the unit to remove any buildup of dust and grime on a regular basis to keep your unit running smoothly.

Also, be sure not to stack any items on or in front of your units venting system. This could cause the compressor to work overtime to maintain temperatures.

Clean Door Gaskets

To ensure that cold air is not escaping, clean any debris or ice build-up from the door gasket to keep an air-tight seal.

Freezer Thermometer

Protect frozen and chilled samples from potential spoilage

Monitor temperatures inside your freezer or refrigerator with the added convenience of sound and flashing LED lights to alert out-of-range temperatures

- Suction cup attaches the probe inside freezer/refrigerator while the thermometer secures to the outside with hook-and loop tape
- Large digital display shows freezer/refrigerator temperature plus ambient temperature
- Switches toggle for freezer/refrigerator monitoring and °F/°C
- Set freezer range: 14.9° to -9.4°F (-9.5° to -23°C);
Set refrigerator range: 35.6° to 46.4°F (2° to 8°C)
- Includes one AAA battery

Item No.	L x W x H in	L x W x H cm
HS28630	4.1 x .79 x 0.9	1.4 x 1.9 x 2.8

Probe Cable measures 0.9 m, Probe measures 5 mm OD x 17 mm L

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